

CARE & MAINTENANCE
DICROMAT FLOW THROUGH PROBE

NOTE: The DiCromat II Salt Analyzer will not provide accurate results if the probe is dirty. It is imperative that it be cleaned. Frequency of cleaning is based on volume of use, as well as type of products. **(We recommend cleaning, at minimum, 1x per week; 1x per shift is preferable.)**

THE FLOW THROUGH PROBE CAN BE CLEANED WITH A SOLUTION OF DAWN DISHWASH DETERGENT AND THE BRUSH SUPPLIED WITH THE UNIT. AFTER CLEANING, THE PROBE SHOULD BE RINSED THOROUGHLY. **ONCE A WEEK, WE RECOMMEND THAT THE PROBE BE CLEANED WITH EITHER AJAX OR COMET CLEANSER.** AGAIN, RINSE THOROUGHLY AFTER CLEANING. REFER TO THE USER'S MANUAL FOR ADDITIONAL INFORMATION ON THE CARE AND MAINTENANCE OF THE FLOW THROUGH PROBE.